

# Dilucup<sup>®</sup> Elegance NaCl (Saline solution)

Ready-to-use dilution cups with 9 mL of sterile buffered diluent.

**Application:** To be used for diluting microbiological samples together with a dedicated shaker, Dilushaker or Dilugent<sup>®</sup> Shaker. By adding a sample of 1 mL to the first cup, a dilution of 1:10 is achieved. 1 mL from this cup is then transferred to the next to get a dilution of 1:100. The procedure is continued until the required dilution is reached.

## Composition per litre of deionized water:

(according to EN ISO 6887-1:2017):

NaCl 8,5 g

pH: 7,0 ± 0,2 at 25°C

Volume per cup: 9 mL ± 2 %

Sterilization: E-Beam

Material of the cups: PET

**Expiry date:** Unopened bags: 30 months after production.

**Storage:** Unopened bags: +2 to +30°C. Opened bags: +2 to +8°C up to 3 months.

**Packaging:** The cups are delivered in blisters of 7 pre-cut rows of 3 cups (40DENACL3) or 6 cups (40DENACL6). Each box contains 672 cups distributed in 2 bags of 336 cups.

**Weight and volume:** 8,1 kg/box - 0,032 m<sup>3</sup>/box

## Article numbers:

40DENACL3: 2 bags x 16 blisters x 21 Dilucups

40DENACL6: 2 bags x 8 blisters x 42 Dilucups

